## **AE-DRY Range Digital With Drying**

Multipurpose vertical sterilizer with final drying

## Gama AE-DRY

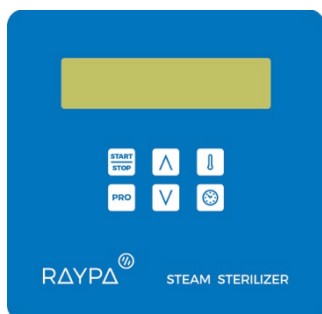
Solids sterilization with multipurpose uses indicated for laboratories and industry. Also suitable for liquids sterilization.

## Operating Principal

AE-DRY vertical range provides a single solution for the multiples sterilization needs on laboratories including instruments, hollowware elements, glassware and other laboratory items that need a final drying to avoid contamination.

Load has to be placed into the vessel's baskets and, after automatic filling with distilled water, device starts to heat up and purge until set temperature is reached. In case of solid objects, discharge is automatically forced during cooling stage to reduce cycle time.

**Digital Microprocessor** – Digital microprocessor for an easy programming and parameters selection for sterilization process.

**Programs** – Microprocessor has 10 programs, from P0 to P9, and the first four are pre-defined and protected.

Program N°	Temperature (°C)	Pressure (bar)	Sterilization time (min.)	Sterilization mode	Drying time (min.)
<b>P0</b>	115	0,6	60	SOLIDS-1	12
<b>P1</b>	121	1,1	30	SOLIDS-1	25
<b>P2</b>	133	2,1	20	SOLIDS-1	30
<b>P3</b>	121	1,1	20	LIQUIDS-1	-

The rest of the programs are user-defined with the following parameters setting:

- Sterilization temperature
- Sterilization time
- Drying time
- Heart probe selection

## Gama AE-DRY

**Water Filling** – Pure water for sterilization can be added manually into the water pan. Automatic filling is also available by connecting the device to pure mains water.

## Sterilization Process

**Pre-vacuum and filling stage** – Integrated vacuum pump for initial pre-vacuum and subsequent automatic water filling.

**Heating Stage** – Fast saturated steam production by means of powerful heating elements.

**Sterilization Control** – A PT-100 temperature probe controls the chamber temperature, with the possibility of installing another flexible PT-100 temperature probe for liquid programs.

**Cooling and drying stage** – In solid program, discharge is automatically forced during cooling stage to reduce the cycle time. Then, integrated heating jacket together with pre-vacuum pulses is activated for a fast and intense drying.

## Gama AE-DRY

**Security** – All safety standards are covered with safety valve, safety thermostat and safety pressure switch, door opening prevention system in case of overpressure, open door sensor and water level detector. EU Directive and Standards:

**2014/35/UE** *Low Voltage*

**2014/30/UE** *Electromagnetic Compatibility*

**2014/68/UE** *Pressure Equipment*

**EN-61010-1**

**EN-61010-2-040**

**EN-61326**

**AD 2000 Merkblatt**

## Communications and Documentation

**RS-232 Port** – For live monitoring of sterilization cycle directly to PC. It is needed the customized software SOFT-S7000W.

**Integrated Printer** – For quality printouts after each cycle. Intervals can be adjustable. Printer has to be requested before ordering.

## Gama AE-DRY

### Technical Data

Referencia Reference Référence	Capacidad l Capacity l Capacité l	Capacidad útil l Usable capacity l Capacité utile l	Ø x H mm	L x D x H mm	Temp. °C	Presión máxima bar Max. pressure bar Pression max. bar	W	Kg	Tiempo min. Time min. Temps min.	Depósito H <sub>2</sub> O l Container H <sub>2</sub> O l Réservoir H <sub>2</sub> O l
<b>AE-28 DRY</b>	33	28	300 x 400	560 x 510 x 1120	100-134	2,1	2000	90	1' - 250'	7
<b>AE-50 DRY</b>	55	50	300 x 700	560 x 510 x 1300	100-134	2,1	3200	110	1' - 250'	7
<b>AE-75 DRY</b>	79	75	400 x 600	730 x 620 x 1190	100-134	2,1	3200	140	1' - 250'	14
<b>AE-110 DRY</b>	115	110	400 x 850	730 x 620 x 1440	100-134	2,1	4500	180	1' - 250'	14
<b>AE-150 DRY</b>	175	150	500 x 750	850 x 760 x 1400	100-134	2,1	6000	265	1' - 250'	24

Electrical connection for equipment of 28, 50 and 75 L: 230V (1P + N), 50/60 Hz.

Electrical connection for 110 and 150 L equipment: 400V (3P + N), 50/60 Hz.

Possibility of modifying the voltage on request.

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#### Intended use

- To lift heavy materials to put into the autoclave.
- Lift with hoist powerful system with adjustable arm.

#### Features

- Easy to use.
- Suitable for vertical autoclaves: 110L and 150L, consult for other models.
- Rotation up to 200 degrees.

#### Safety

- Emergency Stop Button.
- Motor with auto braking system.

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#### SET AUTOMATIC FILLING

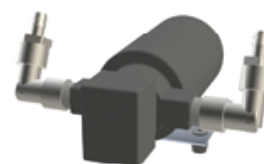
• To be installed in our facilities.

References:  
**KLL:** for B and DRY from 28 up to 150 L  
**KLL-21:** for AH-21B and AH-21L  
**KLL-MP:** for MP

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#### FLEXIBLE "HEART" PROBE PT-100

- Recommended for sterilizing liquids in flasks of 1 litre or bigger.
- The control is done from the microprocessor.
- This accessory should be installed in factory.

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#### THERMAL PAPER PRINTER

- It prints program number, cycle number, temperature, date and hour of the run.
- Error messages are also printed.
- To be installed in our facilities.

References:  
**IT:** LCD Screen  
**IT/TS:** Touch Screen

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Referencia: **ITS**

#### TABLE TOP PRINTER

- It prints program number, cycle number, temperature, date and hour of the run.
- Error messages are also printed.
- It includes RS232 connection.

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\* Serie AES, AE-DRY

#### PC CONNECTION SW7000

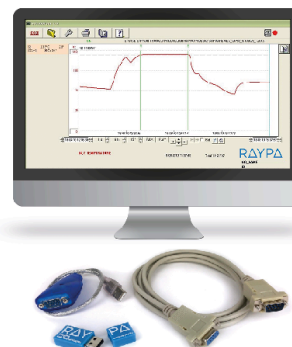
- Communication software between the equipment and the PC.
- Connection for RS-232.
- USB Adapter.
- Not compatible with touch screen circuits.

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\* Todos los modelos / All models / Tous les modèles

#### STERILIZATION CONTROL TAPE

- Chemical indicator for steam. The change of color indicates that the materials have been processed, without being a guarantee of proper sterilization. Class 1.
- To verify that sterilization is effective, additional methods are needed as biological indicators EN ISO 11138.
- Tape roll of 50 m x 19 mm

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