AES Range Semiautomatic Autoclaves

Multipurpose vertical sterilizer

AES Range


Solids and liquids sterilizer with multipurpose uses indicated for laboratories and industry.

Operating Principal

AES vertical range provides a single solution for the multiples sterilization needs on laboratories including liquids, culture media, biological waste, contaminated media, instruments, glassware and other laboratory items.

Load has to be placed into the vessel's baskets and, after manual filling with distilled water, device starts to heat up and purge until set temperature is reached. In case of solid objects, discharge can be manually forced during cooling stage to reduce cycle time.

Digital Microprocessor – Digital microprocessor for an easy programming and parameters selection for sterilization process.

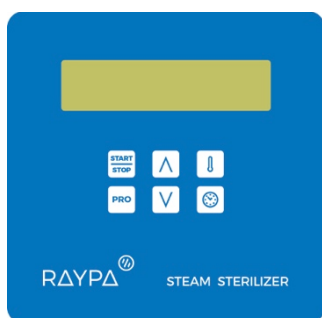Programs – Microprocessor has 10 programs, from P0 to P9, and the first four are pre-defined and protected.

Program N°	Temperature (°C)	Pressure (bar)	Sterilization time (min.)	Sterilization mode
P0	115	0,6	60	SOLIDS-1
P1	121	1,1	30	SOLIDS-1
P2	133	2,1	20	SOLIDS-1
P3	121	1,1	20	LIQUIDS-1

The rest of the programs are user-defined with the following parameters setting:

- Sterilization temperature
- Sterilization time
- Heart probe selection

AES Range

Sterilization Process

Heating Stage – Fast saturated steam production by means of powerful heating elements.

Sterilization Control – A PT-100 temperature probe controls the chamber temperature, with the possibility of installing another flexible PT-100 temperature probe for liquid programs.

Cooling Stage – In solid program, discharge can be manually forced during cooling stage to reduce the cycle time.

Security – All safety standards are covered with safety valve, safety thermostat and safety pressure switch, door opening prevention system in case of overpressure, open door.

EU Directive and Standards:

2014/35/UE *Low Voltage*

2014/30/UE *Electromagnetic Compatibility*

2014/68/UE *Pressure Equipment*

EN-61010-1

EN-61010-2-040

EN-61326

AD 2000 Merkblatt

Communications and Documentation

RS-232 Port – For live monitoring of sterilization cycle directly to PC. It is needed the customized software SW7000.

Integrated Printer integrated – For quality printouts after each cycle. Intervals can be adjustable. Printer has to be requested before ordering.

AES Range

Technical Data

Referencia Reference Référence	Volumen total / útil Total volume / usable Volume total / utile l	Dimensiones útiles cámara Usable chamber dimensions Dimensions utile chambre Ø x H mm	EXTERIOR L x D x H mm	Temp.°C	Presión máxima Max. pressure Pression max. bar	W	Kg
AES-28	33/31	300 x 440	505 x 580 x 1110	100 - 134	2,1	2000	75
AES-50	55/50	300 x 710	505 x 580 x 1290	100 - 134	2,1	3200	90
AES-75	79/75	400 x 600	610 x 700 x 1185	100 - 134	2,1	3200	115
AES-110	115/110	400 x 850	610 x 700 x 1435	100 - 134	2,1	4500	135
AES-150	175/153	500 x 780	750 x 820 x 1400	100 - 134	2,1	6000	230

Electrical connection for equipment of 28, 50 and 75 L: 230V (1P + N), 50/60 Hz.
Electrical connection for 110 and 150 L equipment: 400V (3P + N), 50/60 Hz.

Possibility of modifying the voltage on request.

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Intended use

- To lift heavy materials to put into the autoclave.
- Lift with hoist powerful system with adjustable arm.

Features

- Easy to use.
- Suitable for vertical autoclaves: 110L and 150L, consult for other models.
- Rotation up to 200 degrees.

Safety

- Emergency Stop Button.
- Motor with auto braking system.

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FLEXIBLE "HEART" PROBE PT-100

- Recommended for sterilizing liquids in flasks of 1 litre or bigger.
- The control is done from the microprocessor.
- This accessory should be installed in factory.

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THERMAL PAPER PRINTER

- It prints program number, cycle number, temperature, date and hour of the run.
- Error messages are also printed.
- To be installed in our facilities.

References:
IT: LCD Screen
IT/TS: Touch Screen

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Referencia: **ITS**

TABLE TOP PRINTER

- It prints program number, cycle number, temperature, date and hour of the run.
- Error messages are also printed.
- It includes RS232 connection.

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